



USB Microscope Directions

- Log in to the server.
- Attach microscope to computer using the USB cord
- Open MacIntosh HD
- Open USB Shot Folder and doubleclick the USB Shot icon
- To create a folder to save your images, click on the folder icon
- Click New Folder,  name it, click Create, and click Choose
- Click on Pop-up Menu and choose classroom or student folder
- Turn microscope light on by moving the button to A
- Focus on selected specimen
- Click the button to capture your image. It is saved in your folder

COOL! SO NOW WHAT??

These images can be used in numerous applications - AppleWorks, Inspiration, Kid Pix, Kid Works Deluxe, HyperStudio, etc., etc.

HINTS:

- The stand may be used to help "steady" or stabilize the microscope.
- Use the back light if needed.
- There are four lenses - 1x, 50x, 100x, and 200x. You might want to experiment with the different powers to pick which is best for your project.
- Make sure to set your computer at Thousands of Colors.
- Another way to capture and image, is to click the camera icon  on the screen.